
Repair Instructions

Premium cabinets top panel replacement



To suit: TPR/TLR/TRI/TBW/TUF cabinets 360/520/750/950/1100/1150/1500

Approvals		
Written By:	Reviewed By:	Approved By:
Duncan Smeaton	Stephen Watt	Duncan Smeaton

Safety Reminder


These instructions do not purport to address all potential safety issues, if any, associated with the product's use. It is the responsibility of the user of these instructions to establish appropriate safety and health practices and determine the applicability of regulatory limitations before use.

Before attempting any of the following, perform the actions listed below:

- Turn OFF power to the machine.
- Unplug the machine.

Summary

These instructions are for removing and replacing a top front panel on a TPR/TLR/TRI/TBW/TUF cabinets 360/520/750/950/1100/1150/1500.

Tools Required	
No hand tools required	
Safety Step. Example Shown	
Additional personnel particularly for larger cabinets.	

Kit Contents		
Item	Description	Quantity
N/A	Replacement top panel.	1

Section 1: Preparation

Step 1 - Turn off the power and remove the plug from the outlet.

Step 2 - Place safety step if needed.

Section 2: Remove

Step 1 - Slide the panel up which will free it from the locating blocks on the cabinet.



Step 2 - Pull the panel forward free of the blocks.



Section 3: Installation

Step 1 - Place the panel slightly above its final position matching the rectangular holes with the locating blocks.



Step 2 - Push the panel forward onto the blocks then slide it down to lock into place.



Step 3 - The panel should now be in place and even on all edges.



Section 4: Power ON

Step 1 - Plug the cabinet back in and turn on the power.

Section 5: Support and Contact

Repair and Support is available over the telephone Monday through Thursday from 8:30am to 4pm and Friday 8:30am to 2pm. Please contact service@thermoline.com.au for email technical support.

You can also visit our website at www.thermoline.com.au for access to additional useful troubleshooting guides, operating manuals, and technical information.

Thermoline Scientific Equipment
10-12 Ross Place
Wetherill Park NSW 2164
Ph: 02 9604 3911
Email: hello@thermoline.com.au
Web: www.thermoline.com.au

