

---

## Repair Instructions

---

---

To replace the door gasket in a TPR145/200 Pharmacy Fridge

---



To suit: TPR/TLR 145 and 200

Approvals		
Written By:	Reviewed By:	Approved By:
Darren Assey	Duncan Smeaton	Duncan Smeaton

## Safety Reminder

These instructions do not purport to address all potential safety issues, if any, associated with the product's use. It is the responsibility of the user of these instructions to establish appropriate safety and health practices and determine the applicability of regulatory limitations before use.

Before attempting any of the following, perform the actions listed below:

- Turn OFF power to the machine.
- Unplug the machine.

## Summary

Instructions on how to remove a old door gasket and install a new one.

Tools Required	
None required	

Kit Contents		
Item	Description	Quantity
WRA21255	Door Gasket to suit TPR-145	1
or		
WRA29931	Door Gasket to Suite TPR-200	1

For the purpose of this guide, the fridge shown in the pictures has is a TPR/TLR-200. However, the process for removing and installing the door seal for a TPR/TLR-145 fridge is exactly the same.

## Section 1: Preparation

Empty the fridge of any stock (maintaining any cold chain requirements) and turn off the refrigerator and remove the plug from the power point.

## Section 2: Remove the old door seal



Open the fridge door and from the top corner, pinch the door seal and begin to peel it away.



Slowly work your way around the door, removing the door seal from the groove until it is completely detached.

### Section 3: Installation



Start from the top corner of the fridge, firstly orientate the new seal to the door and then take the corner of the seal and push it into the groove of the door frame. Make sure the ridged side of the door seal is the side that is inserted into the groove.



Push the seal right along each edge of the door until it is fully inserted. Check to see if it is seated correctly.

## **Section 4: Power On**

Plug the fridge back in and turn on the power. Wait for the Fridge to get to temperature before replacing the stock.

## **Section 5: Support and Contact**

Repair and Support is available over the telephone Monday through Thursday from 8:30am to 4pm and Friday 8:30am to 2pm. Please contact [service@thermoline.com.au](mailto:service@thermoline.com.au) for email technical support. You can also visit our website at [www.thermoline.com.au](http://www.thermoline.com.au) for access to additional useful troubleshooting guides, operating manuals, and technical information.

Thermoline Scientific Equipment  
10-12 Ross Place  
Wetherill Park NSW 2164  
Ph: 02 9604 3911  
Email: [hello@thermoline.com.au](mailto:hello@thermoline.com.au)  
Web: [www.thermoline.com.au](http://www.thermoline.com.au)

